

Baths, Uncirculated (Ambient +5 °C to +100°C)

AUSTRALIAN MADE

| ORDER CODE | DIM. INT USEABLE LxWxD cm | DIM. EXT LxWxD cm | LITRES CAPACITY TOP - 4cm | ELECT. 240V 50Hz | APPROX WEIGHT Kg | OPTIONAL POLYCARB. LID | OPTIONAL RESET THERMOSTAT |
|------------|---------------------------|-------------------|---------------------------|------------------|------------------|------------------------|---------------------------|
| TWB-10 | 23x30x17 | 35x32x22 | 10 | 500W | 8 | PCL-10 | MRST-10 |
| TWB-22 | 50x30x17 | 61x32x22 | 22 | 900W | 11 | PCL-22 | MRST-22 |
| TWB-40 | 58x48x17 | 72x51x22 | 40 | 2000W | 15 | N/A | MRST-40 |

FEATURES

- Insulated Uncirculated Bath standard sizes 10, 22, and 40 litre capacity
- Available in Thermostat Control “T” or Digital Control “D” with Display
- Seamless Stainless Steel tank (eliminates bug growth in corners)
- Radial corners (quick and easy to clean or disinfect)
- Carry handles fitted as standard
- Perforated removable floor
- Custom baths can be ordered to suit your application
- Temperature range Ambient +5°C to +100°C



TWB-10D

Caution:

The **Optional** Manual Reset Safety Thermostat (MRST) is a fail safe device factory set to cut out the heating if the bath runs dry. This is particularly important if the bath is operated whilst unattended for long periods, or at elevated temperatures.

| HEAT UP TIMES (min) | AMB. TO 37°C | AMB. TO 50°C |
|---------------------|--------------|--------------|
| TWB-10 | 30 | 45 |
| TWB-22 | 40 | 60 |
| TWB-40 | 50 | 90 |

Temperature Range:

ANALOGUE: +30°C to +100°C
 DIGITAL: Ambient +5°C to +100°C
 CONTROL STABILITY: ±0.25°C

Water Bath Polycarbonate Floating Ball Blanket.

Prevents evaporation and helps maintain constant bath temperature.

| | |
|---------------|---|
| TPB 10 | Bag of 200 balls to suit 10 Litre bath. |
| TPB 22 | Bag of 400 balls to suit 22 Litre bath. |
| TPB 40 | Bag of 750 balls to suit 40 Litre bath. |



Optional Accessories

- ✓ Polycarbonate gable lid for the TWB-10 and TWB-22 Series
- ✓ Stainless Steel gable lid for the TWB-40 series
- ✓ Polycarbonate Floating Balls [replaces above lid] and prevents evaporation
- ✓ Safety Devices optional Manual Reset Safety Thermostat Device (MRST) prevents damage to bath and bench should the bath water level fall to a dangerous level during long periods of unattended operation.